


Rapid Separation of Basic Drug Compounds Under Alkaline pH Conditions on pH-Stable, Hamilton PRP™-C18

Mobile phase pH is a powerful tool in methods development, particularly for separation of neutral forms of amines or other organic bases under alkaline conditions. In this study a generic, 5-min linear gradient was used to separate six basic drug compounds on a short (50 mm) PRP-C18 column.

More than 70% of all pharmaceutical drug compounds are cationic solutes that carry a formal positive charge below pH 7. Separation of these and other organic bases has historically been problematic. Ionization has a dominating effect in reversed phase chromatography that tends to dictate retention. Consequently, the elution window for a sample of ionized amines is narrow. The task is further complicated by secondary interactions that occur between positively charged solutes and residual silanols on the column stationary phase. These secondary mechanisms of retention are the principle source for anomalous chromatographic activity, such as poor peak shape, shifts in retention times and loss of efficiency that progressively worsen over the life of the column.

The PRP-C18 is a new column designed for high-efficiency reversed phase separations under any mobile phase conditions. The stationary phase for the PRP-C18 is devoid of free silanols, does not strip, bleed, or dissolve at any pH, and therefore can be expected to perform reliably and reproducibly throughout the extended life of the column, regardless of mobile phase conditions. Use of alkaline mobile phase (pH > 11) permits separation of basic solutes in their neutral forms. This broadens the window for elution, whereby subtle structural nuances among chemically similar compounds can be exploited to effect resolution.

Although some recent C18 columns boast stability in alkaline pH, all silica-based supports experience measurable degradation at pH > 6, where column life is still considerably shorter than if used under more favorable conditions. On the other hand, the PRP-C18 stands up to prolonged exposure to concentrations as high as 1 M NaOH and H₂SO₄, with no measurable decrease in performance.

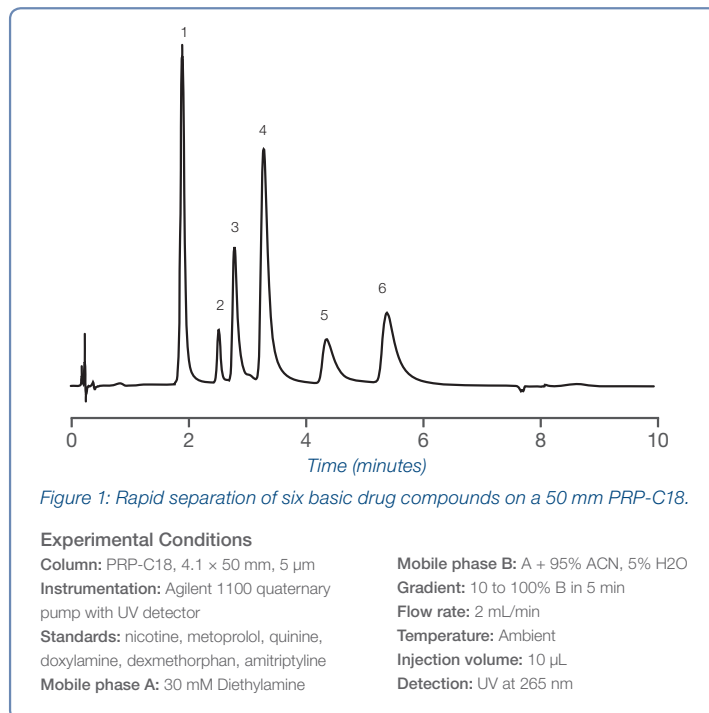

Results and Conclusion

In modern drug discovery science where analytical HPLC can be a bottleneck, the trend is to streamline production through the use of shorter columns with smaller particles operated at elevated flow rates. The flexibility to employ a high pH mobile phase is another valuable tool that permits separation of basic solutes in their neutral forms. Often times this greatly simplifies the process of methods development. In this study, separation of a set of structurally diverse pharmaceutical compounds is achieved on a short (50 mm) PRP-C18 column using a generic 5 min linear gradient.

Authors:

Derek Jensen and Mark Carrier, Hamilton Company

© 2013 Hamilton Company. All rights reserved.
All trademarks are owned and/or registered by Hamilton Company in the U.S. and/or other countries. Lit. No. L80091

HAMILTON®

Web: www.hamiltoncompany.com

USA: 800-648-5950

Europe: +41-81-660-60-60

Hamilton Americas & Pacific Rim
Hamilton Company Inc.
4970 Energy Way
Reno, Nevada 89502 USA
Tel: +1-775-858-3000
Fax: +1-775-856-7259
sales@hamiltoncompany.com

Hamilton Europe, Asia, & Africa
Hamilton Bonaduz A.G.
Via Crusch 8
CH-7402 Bonaduz, Switzerland
Tel: +41-81-660-60-60
Fax: +41-81-660-60-70
contact@hamilton.ch

To find a representative in your area, please visit hamiltoncompany.com/contacts.